



704

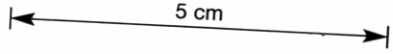

## Nomogram for Specific Gravity & Temperature Correction in Cyclosizer

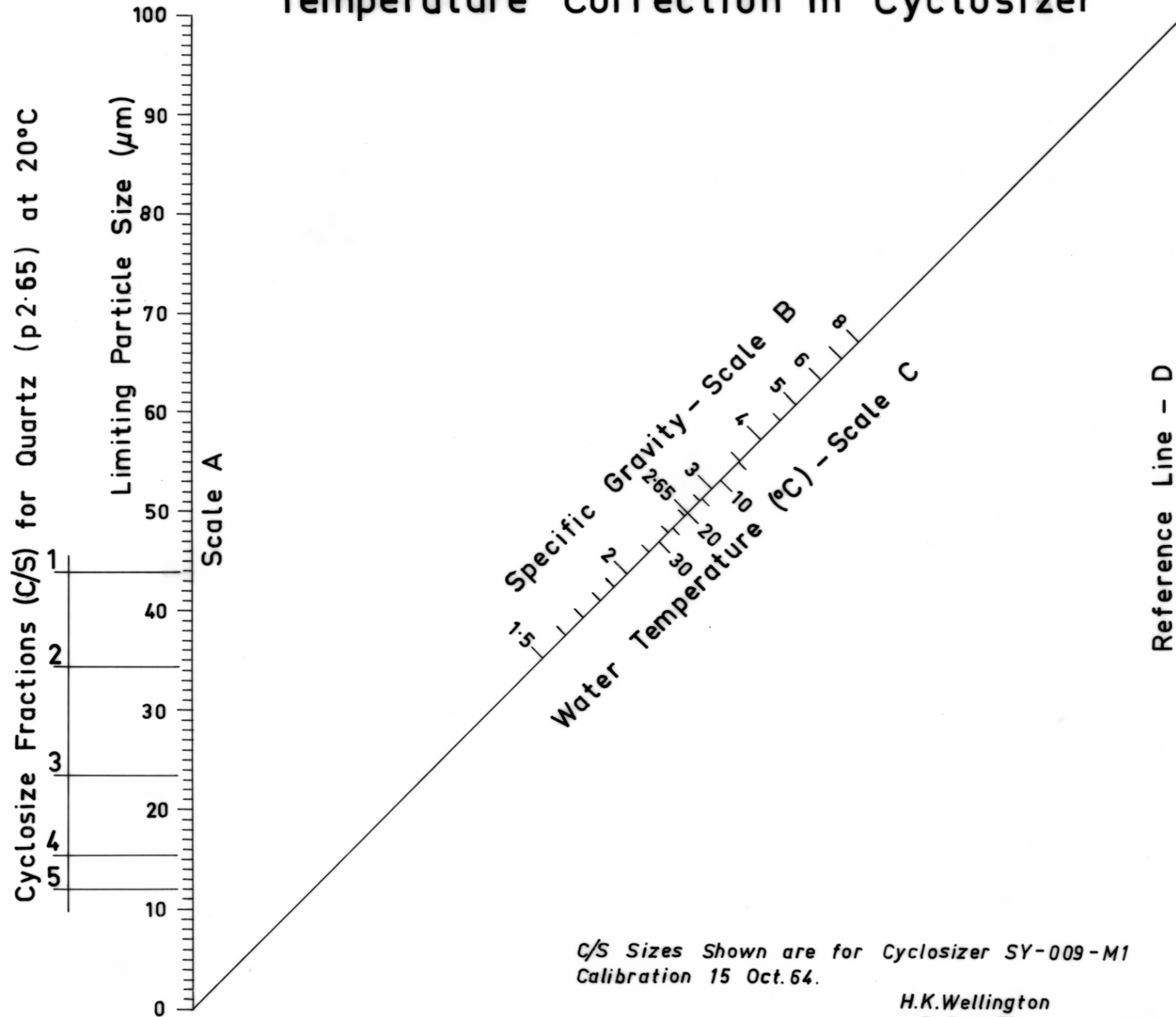

*How to use :*

1. Connect C/S Fraction Size on Scale A with specific gravity of Scale B and note the intercept on the Reference Line D
2. Connect intercept on Reference Line D with water temperature on Scale C and read the corrected size on Scale A.

*Example :*  
 C/S 2 for SG4 @ 10°C  
 is 29 $\mu\text{m}$

*C/S Sizes Shown are for Cyclosizer SY-009-M1  
 Calibration 15 Oct. 64.*

H.K. Wellington  
 7 Sept. 72